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Instruction for Use TPP Serological Pipette

TPP's sterile serological pipettes are for dispensing aqueous fluids such as cell culture media.

The color-coding on the pipette and dispenser allows for easy identification.

The negative graduation and reservoir provide more working volume than conventional serological pipettes.

The specially designed dispenser with perforated peel-off tab serves as a dispenser for easy collection directly at the bench.

The TPP serological pipette is for single use only.

Note:

Follow the national regulations for handling biological materials and wear the appropriate protective clothing.

During the work process, be sure to observe the rules of aseptic technique.

Do not use your mouth to pipette; always use the pipetting aid to fill and to dispense.

Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use.*

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Advice and Recommendations:

- *If pipetting in a sterile environment and the pipette touches a non-sterile surface, discard the pipette and replace it with a sterile one.*
- *Risk of contamination with multiple use and/or solution change of pipette: Always replace the pipette when changing solutions or cell suspensions to avoid cross-contamination.*

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Chemical Resistance Ink

+ = Good chemical resistance Substance does not cause damage	± = Good to limited chemical resistance Long-time and/or continuous exposure to the substance causes minor damages	-- = Poor chemical resistance Not suitable for exposure to the substance, damage may occur (deformation, discoloration, cracking, dissolution).
--	--	---

Solution	+ / ± / -	Solution	+ / ± / -
Acetic acid	±	Formaldehyde	±
Acetone	-	Glutaraldehyde	-
Acetonitrile	-	Glycerol	+
Ammonia solution	±	Isoamyl alcohol	±
Caustic soda	-	Isopropanol	±
Chloroform	-	Methanol	+
Dimethyl sulfoxide	+	Nitric acid	-
DMEM	+	Sulfuric acid	-
Ethanol	±	Triton X-100	+
Fetal bovine serum	+	Water	+

Verpackungseinheit 3-fach verpackte TPP Serologische Pipetten / Packaging unit for triple bagged TPP Serological Pipettes

Nr. No.	Karton Case	Produktbeschreibung 3-fach verpackt Product Description Triple Bagged	Stk. / Beutel Qty. / Bag	Stk. / Karton Qty. / Case
394525	1/4	Serologische Pipette 25 ml mit Reservoir <i>Serological pipette 25 ml with reservoir</i>	1 bag of 3	36
394550	1/4	Serologische Pipette 50 ml mit Reservoir <i>Serological pipette 50 ml with reservoir</i>	1 bag of 3	36

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Additional:

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from www.tpp.ch.